



Xylene Isomers (99% p-Xylene)

Column: **CHFFAP - 50m, 0.32mm, 1.00µm**

Catalog Code: CHFFAP-5032100

Conditions:

Injection: Split 250°C, 0.5µL, 350mL/min Split Flow.

Detector: FID 250°C.

Oven Temp.: 65°C, 2°C/min, 150°C.

Carrier Gas: Helium, 150 kPa.

Peaks:

1. Ethyl Benzene
2. p-Xylene (99%)
3. m-Xylene
4. o-Xylene

For more information, please contact:

info@chrom4.com | www.chrom4.com